



ULTRA-SONIC CLEANER #2

TARTAR AND STAIN REMOVER

Ultra-Sonic Cleaner #2 is specially formulated to remove mineral scale, surface deposits, and stubborn stains from a variety of hard, non-porous items. Effective in ultrasonic cleaning systems for laboratory, industrial, maintenance, and hobby applications.

—Removes mineral deposits and stains from hard surfaces

—Fast working

Indication for use: Removes mineral scale and stains from hard, non-porous items. Not for medical use.

Standard Size Available: 1 Gallon (3.8 Liters)

Directions: Place item in covered glass beaker of undiluted Ultra-Sonic Cleaner #2. Immerse the covered glass beaker in Ultra-Sonic Cleaner #1 Solution. USC-1 acts as a conductor between the tank and the beaker. DO NOT USE ON ALUMINUM. Cleans in 5-10 minutes. Rinse items with water after removing them from Ultra-Sonic Cleaning Solution.

Item #: 3590040



RELATED PRODUCTS

Ultra-Sonic Cleaner #1: General Purpose Concentrate - 3570040

Ultra-Sonic Cleaner #5: Plaster and Stone Remover - 3610040

Manufactured by Garreco, LLC
430 Hiram Road, Heber Springs, AR 72543, USA
Toll Free: (800) 334-1443
garreco.com

Garreco, LLC is an ISO 13485 Facility
Garreco is a trademark of Garreco, LLC, Heber Springs, AR. ©2025 All rights reserved.